



Vesicle Flow Cytometry Analysis Kit

FOR VESICLE COUNTING AND SIZING

Instrument Setup:

Beckman Coulter CytoFlexS™

Instrument Setup - Beckman Coulter CytoFlex™

Prior to running the vFC Assay, configure the CytoFlex with the appropriate filters and create an Acquisition Template. Typically, this will only need to be performed once, and the saved Template and Experiments used for future experiments.

Filter configuration. Configure the CytoFlex™ for Violet Side Scatter (VSSC) detection. If using a core instrument you will need permission from the core manager. A detailed description of how to perform this operation can be found in the CytoFlex Operation Manual. Briefly, you will access the Violet laser optical filters inside the instrument, place the Violet 405nm filter in position 2, the Violet 450nm filter in position 3, and an unused filter in position 1.

Acquisition Template. The CytoFlex acquisition settings are stored in a .xitm file, which contains data acquisition setting as well as the file and group naming rules applied to file names and Sample IDs. Frequently used plate layouts can be saved and used as templates for future experiments.


To configure an initial Acquisition Template (.xitm file):

1. Configure detectors
2. Define plate and sample wells
3. Set Acquisition conditions
4. Adjust Trigger channel threshold

For subsequent runs, recall the vFC Assay Template and:


1. Define a new Plate and Sample wells
2. Set new File and Group Names

3. Set Acquisition conditions.

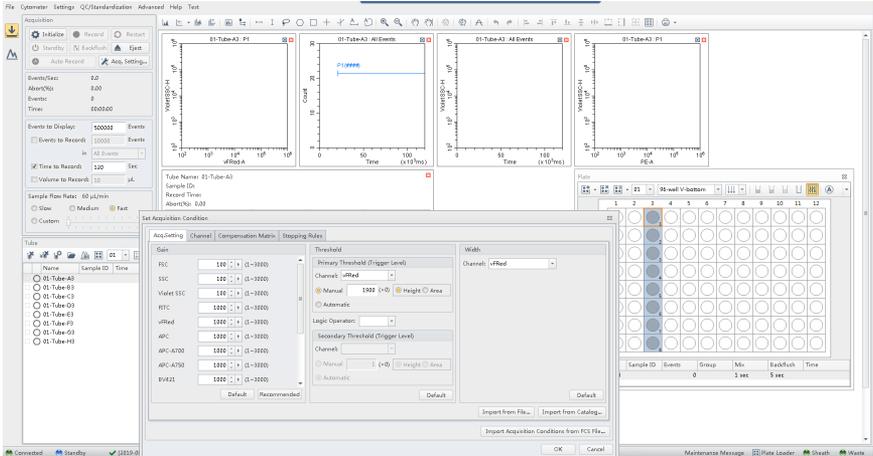

On the *Acq.Setting* tab, set *Gains* for all fluorescence channels at 100, 488 nm scatter (*FSC* and *SSC*) at 100, 405 nm scatter (*VSSC*) at 100.

Set the *Primary Threshold* to the vFRed channel and select the Height (-H) parameter and set the *Manual* threshold to 1900. Set the *Width* channel to vFRed (690/50).

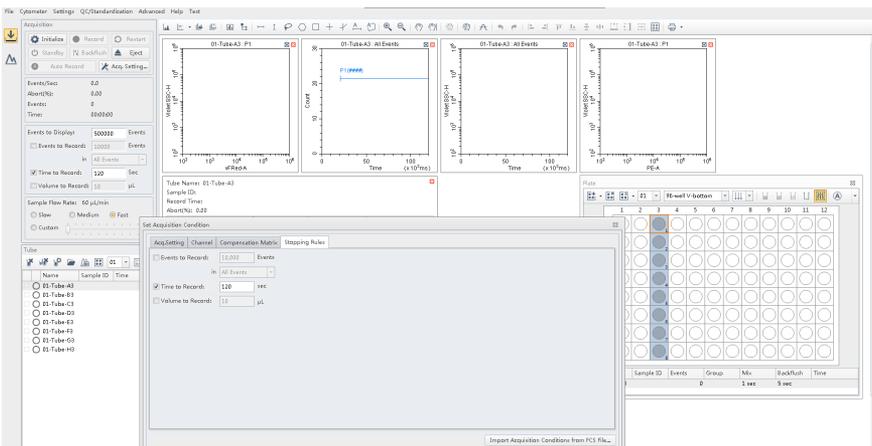

On the *Stopping Rules* tab, check *Time to Record* and enter 120 seconds. Uncheck *Events to Record*. At High Sample Speed (~60 uL/min) on the CytoFlex, this will collect data from 120 ul of sample.

Note: Typically, data from the first 20 ul (20 seconds) are discarded, due to high background that frequently occurs at the beginning of a well, so that data from 100 uL is analyzed for each sample. This gating will happen automatically using the vFC™ Data Analysis Layout.

In the *Acquisition* dialogue, set *Events to Display* to 500,000.

4. **Adjust Trigger Channel Threshold.** Load a well of VFC Staining Buffer and allow system to run for 15 seconds on high flow rate (60ul/min). Set the instrument to trigger from the blue laser-excited, 690/50-H (vFRed) fluorescence channel at 1900, then adjust the manual threshold until you get 10-15 events/second in a Buffer Only sample. Apply this threshold to all wells and save template file (.xitm) for future use. In general, once optimized the threshold does not need to be changed from day to day.

Save Template as vFC Assay.

5. Set File and Group names.

Table from 'Set Naming Rules' dialog:

Name	Sample ID	Events	Group	Mx	BackFash	Time
01-20190911 protoco...	Lipo100 size stand...	37915	Lipo100 size stand...	3 sec	3 sec	2019-09-11 13:...

Highlight wells and define Names as yearmonthdate (eg. 20190704) and Set Group names as defined in “staining wells” on the relevant Protocol Plate Map.

Table from 'Set Group' dialog:

Name	Sample ID	Events	Group	Mx	BackFash	Time
01-20190911 protoco...	buffer only	3813	buffer only	3 sec	3 sec	2019-09-11 13:...

6. Auto Record data. You are now ready to run samples. Click Auto Record to run the selected Sample wells.

Notes